

Imagen® 50-1500bp DNA Ladder

Cat. No DL001

Catalog Number	Size	Concentration
DL001	100 reactions	56 µg / 500 µl

Description


Imagen® 50-1500bp DNA Ladder is combined with PCR products and lots of proprietary plasmids digested with appropriate restriction enzymes to yield different fragments. This product is suitable for use as molecular weight standards ranging from 50-1,500 base pairs for agarose gel electrophoresis. The 200, 500 and 1200 base pair bands have increased intensity to serve as reference points. The approximate mass of DNA in each band can estimate by the intensity of similar size band of ladder (0.56 µg with 5µl loading).

Source

PCR products and double-stranded DNA digested with appropriate restriction enzymes, are phenol extracted and equilibrated to 10 mM Tris-HCl (pH 8.0) and 10 mM EDTA.

Storage

- ✓ Stable for up to 6 months at 4°C.
- ✓ Stable for up to 24 months at -20°C.

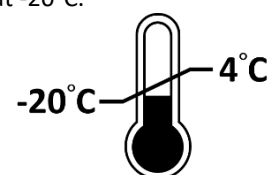


Reaction Setup

5 µl / well, containing orange G as tracking dye.

Application

- ✓ Gel electrophoresis
- ✓ Range: 50-1,500 bp

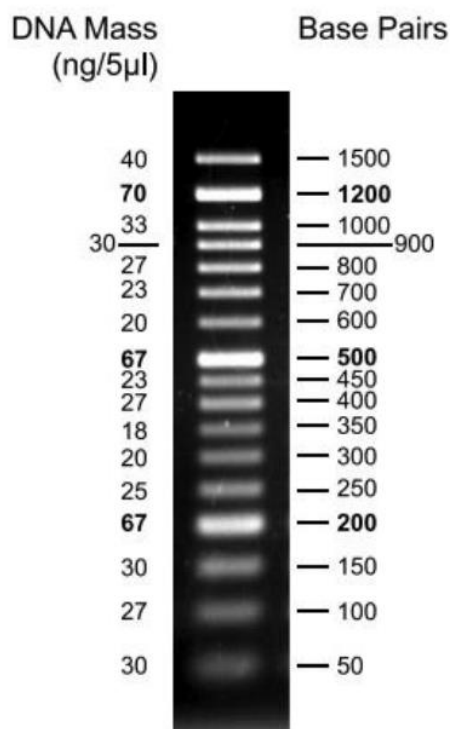

2% TAE agarose gel